



GUARDIAN™ SERIES
Hotplate Stirrers

Think Safer. Think SMART.

Modernize your lab with the
Guardian Series Hotplate Stirrers
designed with outstanding safety and
superior performance in mind.



Ingenuously Practical

Think Safer. Think SMART.

Modernize Your Lab with the Guardian™ Series Hotplate Stirrers

Featuring OHAUS SMART technology, the Guardian Series protects the lab with advanced safety features designed to reduce the chance of accidents. Designed with superior heating and stirring performance, the Guardian 7000, 5000, 3000, and 2000 Series Hotplate Stirrers combine durability and affordability to reduce the high costs of doing business in science labs.

Modernize your lab with OHAUS SMART technology including ***SmartPresence™**, ***SmartLink™**, ***SmartRate™**, ***SmartHeat™**, **SafetyHeat™**, and ***SmartHousing™**. The bright LCD screen displays temperature, speed and *time, and a green indicator lights when the unit is heating and stirring.

Contents


Guardian Series	1
Guardian 7000	2
Guardian 5000	4
Guardian 3000	6
Guardian 2000	8
Frequently Asked Questions	10
Comparison Chart	12
Accessories	13


(*Smart Features, SafetyHeat and Timer are model specific. See page 12 comparison chart.)


GUARDIAN™ 7000


Durable Hotplate Stirrers with Superior Safety Features and Intelligent Performance


OHAUS Guardian 7000 Hotplate Stirrers are designed with outstanding safety features and superior heating and mixing performance. The SmartHousing is chemical resistant with an innovative, easy-to-clean design that channels spills away from internal components and the angled glass control panel. The bright LCD display includes intuitive icons, temperature, speed and time settings - all designed to monitor performance from across the lab.



SmartPresence


SmartLink


SmartHeat


SmartHousing


SafetyHeat


SmartRate



Model e-G71HS07C

Unique Features



SmartPresence™ and SmartLink™ Technologies Ensure Supervised Operation

Protect your lab with patented SmartPresence and SmartLink technologies. These exclusive features automatically shut off the heater if no one is detected after a user set “time out” period has passed or the *Bluetooth® link between a paired mobile device is broken.

SafetyHeat™ System Protects the Lab

An industry-leading early detection system that uses two independent safety controls to continuously monitor the electronics and shut off heating before an overtemperature condition occurs.

Powerful, Consistent Stirring Performance

Precise stirring from 60 to 1600 rpm is made possible by the design of a powerful motor, strong magnet and software controlled ramp rate, ensuring secure magnetic coupling for viscous applications.

Customized Performance for Precise Applications

SmartHeat™ allows you to set the maximum temperature of the unit and SmartRate™ enables fast or slow temperature and speed ramp rates. The programming feature allows storage of 5 multi-step programs.

Model No.	e-G71HS07C	e-G71HS10C	e-G71HSRDM
Top Plate Size	7 x 7 in / 17.8 x 17.8 cm (Square)	10 x 10 in / 25.4 x 25.4 cm (Square)	5.3 in diameter / 13.5 cm (Round)
Capacity	15 L	18 L	20 L
Function	Heating and Stirring		
Top Plate Material	Ceramic		Aluminum
Speed Range	60 rpm to 1600 rpm		
Temperature Range	Ambient + 5 °C – 500 °C		Ambient + 5 °C – 380 °C

*The Bluetooth word mark and logos are registered trademarks owned by Bluetooth SIG, Inc. and any use of such marks by OHAUS is under license.

GUARDIAN™ 5000

Expertly Engineered Hotplate Stirrers with Premium Safety and Versatility Features for Everyday Use

The Next Generation Guardian 5000 Hotplate Stirrers are engineered for safety and versatility across research, academic and industrial settings. They offer powerful heating and mixing, with premium hardware and software to support diverse users safely and reliably. Heating models reach up to 550 °C and feature 3 temperature control modes for quick heating or precise heating effectively reduce temperature overshoot. Select from 5 models with 3 plate sizes.



Model e-G52HSRDA



Unique Features



Safety Features to Protect the Lab

Our signature SafetyHeat™ and SmartHeat™ features combined with the new Run Dry Protection system plus the Time Run mode and the optional Splash Shield provide complete safety.

The Ideal Choice for Various Applications

The evolved SmartRate™ system supports heating & stirring from fast and intensive to soft and precise, with a range of accessories to empower a variety of applications.

Easy to Read LCD Display With Everything You Need to Know

An intuitive user interface featuring a large, bright LCD screen with white characters & knob button(s), allows for user friendly operation. RS232 port included for remote control or data logging.

Model No.	e-G52HSRDA	e-G52HS10C	e-G52HS07C	e-G52HP07C	e-G52ST07C
Top Plate Size	Ø 5.3 in / 13.5 cm	10 x 10 in / 26 x 26 cm	7 x 7 in / 18 x 18 cm		
Capacity	20 L	22 L	20 L	18 L	20 L
Function	Heating and Stirring			Heating	Stirring
Top Plate Material	Ceramic Coated Aluminum	Ceramic			
Speed Range	50 - 1800 rpm			--	50 - 1800 rpm
Temperature Range	Ambient + 5 °C – 360 °C	Ambient + 5 °C – 500 °C	Ambient + 5 °C – 550 °C		--

GUARDIAN™ 3000

Affordable Digital Hotplate Stirrers
Designed for Safety, Powerful
Stirring and Fast Heating

The OHAUS Guardian 3000 Hotplate Stirrers carry the everyday load from classroom to lab.

Protect your lab and classroom with the affordable Guardian 3000 Series Hotplate Stirrers that deliver dependable performance with fast heating, and reliable stirring in everyday applications. The design ensures temperature control of the top plate or your sample when using the optional temperature probe and has a Single Point Calibration feature. All Guardian 3000 Series models offer a bright LCD display that features easy-to-read temperature and speed settings, and intuitive icons.



Model e-G31HSRDS



Unique Features



Never Compromise on Safety

Industry-leading safety at affordable prices. Guardian 3000 units offer safety, controlled heating and powerful stirring for general laboratory use – all at a competitive price point.

SafetyHeat™ System Protects the Lab

SafetyHeat protects the lab. An industry-leading early detection system uses two independent safety controls to monitor electronics and stop heating before overheating occurs.

Reliable and Durable for Every Lab

Designed to last - the rugged metal housing has a spill-resistant design for long life in the lab under harsh conditions. The smooth, chemical-resistant ceramic top makes for easy clean-ups.

Designed to be Safe

Consistently safe to use at high temperature settings with a spill-resistant design that features a rugged metal housing that delivers durability in the lab under harsh conditions. The chemical-resistant ceramic top plate is designed for easy clean-up.

Model No.	e-G31HS04C	e-G31HS07C	e-G31HSRDS
Top Plate Size	4 x 4 in / 10.2 x 10.2 cm (Square)	7 x 7 in / 17.8 x 17.8 cm (Square)	5.3 in diameter / 13.5 cm diameter (Round)
Capacity	15 L		
Function	Heating and Stirring		
Top Plate Material	Ceramic		Ceramic Coated Stainless Steel
Speed Range	80 rpm to 1600 rpm		
Temperature Range	Ambient + 5°C – 500°C		Ambient + 5°C – 380°C

GUARDIAN™ 2000


Rugged and Reliable Analog
Hotplates and Stirrers for General
Purpose Heating and Stirring

Dependable OHAUS Guardian 2000 Hotplates and Stirrers are ideal for education and other lab environments that require affordable, everyday fast heating and reliable stirring up to 15L. All three sizes offer simple knob control to easily adjust the heating and stirring functions. The durable housing includes a built-in support rod holder which accepts the optional support rod. Stirring models include PTFE-coated stir bars.



Model e-G21HS07C

Unique Features

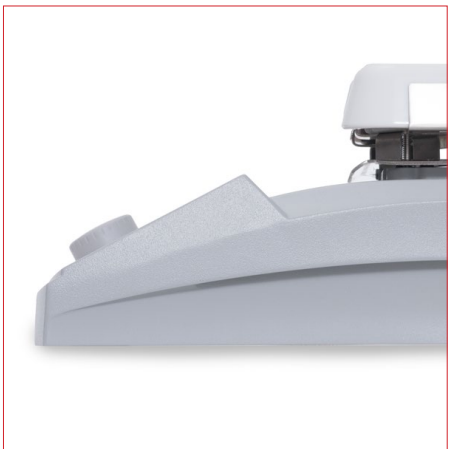

Consistent, Dependable Stirring and Heating

Powerful stirring and fast heating at an economical price. Safely heat up to 500°C and stir up to 15L with an easy-to-use knob control for simple and reliable everyday heating and stirring in the lab.

Designed for Reliability and Durability in Every Lab

Designed to last - the rugged metal housing has a spill-resistant design for long life in the lab, under harsh conditions. The smooth, chemical-resistant ceramic top plate is perfect for easy clean up.

Designed for User Safety

Safety is of utmost importance - the prominent hot top indicator light illuminates when the heater is above 40°C, keeping the end user safe during all heating applications.

Choose From 9 Models to Suit Your Unique Application

The Guardian 2000 Series is available in three top plate sizes: 7x7 inch (17.8 x 17.8 cm), 4x4 inch (10.2 x 10.2 cm) and round top (5.3 in / 13.5 cm diameter). Each size features three models - dual function hotplate-stirrer, and single function hotplate or stirrer.

Model No.	e-G21ST04R	e-G21HP04C	e-G21HS04C	e-G21ST07R	e-G21HP07C	e-G21HS07C	e-G21STRDS	e-G21HPRDS	e-G21HSRDS
Top Plate Size	4 x 4 in / 10.2 x 10.2 cm (Square)			7 x 7 in / 17.8 x 17.8 cm (Square)			5.3 in diameter / 13.5 cm diameter (Round)		
Capacity	15 L								
Function	Stirring	Heating	Heating and Stirring	Stirring	Heating	Heating and Stirring	Stirring	Heating	Heating and Stirring
Top Plate Material	Resin	Ceramic		Resin	Ceramic		Ceramic Coated Stainless Steel		
Speed Range	200 - 1600 rpm	–	200 - 1600 rpm	200 - 1600 rpm	–	200 - 1600 rpm	200 - 1600 rpm	–	200 - 1600 rpm
Temperature Range	–	70 °C - 500 °C		–	70 °C - 500 °C		–	50 °C - 380 °C	

GUARDIAN™ SERIES

Frequently Asked Questions

What differentiates the Guardian 7000 Hotplate Stirrers?

SmartLink™ – provides long-range user detection for improving safety when using the heating function.

SmartPresence™ – detects when a user is nearby and safely turns off the heating function when no one is detected.

SmartHousing™ – ensures that every Guardian hotplate stirrer is consistently safe to use at high temperature settings by remaining cool to the touch. The innovative design is easy to clean and channels spills away from the internal components, as well as being constructed from a chemical resistant material for long-term use in the lab.

How can the Guardian 7000/5000 safeguard sensitive applications?

For precise sample control with the Guardian 7000/5000, use the included temperature probe and enable **SmartRate™** to select fast or slow temperature and speed ramp rates.

SmartHeat™ allows lab users to limit the maximum temperature of the hotplate stirrer preventing overheating of sensitive samples keeping the user and samples safe.

What type of Guardian top plate material best suits the lab? Ceramic, Ceramic Coated Steel, Ceramic Coated Aluminum, Aluminum, or Resin

Ceramic, ceramic-coated steel, ceramic-coated aluminum, aluminum and resin top plates are all suitable for lab use and each material has its benefits.

Ceramic top plates are chemical-resistant and safer if working with corrosive chemicals that may splash onto the plate surface. They can withstand higher temperatures up to 500 °C, and can be wiped clean.

The white surface of a ceramic, ceramic-coated steel or ceramic-coated aluminum top plate is preferable for quantitative chemical analysis or other work where clear visibility of the sample color is required.

The edges of a ceramic top plate may not be as hot as the center. This may not be suitable if uniform heating over the entire plate surface is needed for the application.

Aluminum top plates offer more uniform temperature over the entire heating surface, and will not crack or chip, and will withstand accidental handling. Ceramic-coated steel and ceramic-coated aluminum top plate materials also provide more uniform temperature across the heating surface with a white surface to easily view the sample. These materials have a lower maximum temperature than the full ceramic top plates.

The durable, chemical-resistant resin top plate on stir only models provides a bright white surface for clear visibility. This long-lasting resin material will not chip or crack if mishandled.

To avoid thermal shock, glassware should be used on ceramic top plates and heating of metallic vessels should be avoided. Ceramic top plate units are also not suitable for sandbaths or any vessel that may reflect heat back onto the surface which may cause thermal shock.

What are the benefits of using the Guardian Series Hotplate Stirrers with round top plates?

The Guardian round top models are ideal for biochemistry and chemistry labs that carry out complex heating and stirring applications and increase productivity by utilizing multiple positions to perform experiments in parallel. Round top models save valuable bench space and accept a wide range of aluminum block accessories that can accommodate standard vials and test tubes.

Which safety features are built into the Guardian Series for use in academic settings?

All Guardian heating models feature a prominent 13 mm Hot Top Indicator Light to alert the end user that the top plate surface is hot and not safe to the touch. This hot top light stays on even when the power switch is turned off. This important feature will help prevent accidents in all settings and especially academic labs where students are using hotplates and hotplate stirrers.

The Guardian 2000 models feature simple overheat protection where the unit will turn off the heating function if the temperature exceeds the upper limit.





Guardian 3000, 5000 and 7000 Series heating models feature **SafetyHeat**; an industry-leading early detection system that uses two independent safety controls to monitor the electronics and stop heating before an overtemperature condition occurs. This important feature protects the sample, the user, and the lab by preventing accidents that occur due to overheating samples.

The digital control Guardian 3000, 5000 and 7000 models are designed so the user must press the knob to start the heating and stirring function. This knob press is also needed to change a setting. This secondary step provides an additional layer of safety by making the user acknowledge the setting. The analog control Guardian 2000 are simple knob control and easy to use where turning the knob starts the function or adjusts the current setting.

All Guardian Series models have chemical-resistant housings, and feature an easy-to-clean design that channels spills away from internal components, and the front panel which protects the unit. The Guardian 7000 Series models are constructed with the **SmartHousing™** that remains cool to the touch at all temperature settings. The Guardian 2000/3000/5000 Series feature a rugged metal housing that will withstand harsh conditions in the lab.

GUARDIAN SERIES

Comparison Chart

	Guardian 7000	Guardian 5000	Guardian 3000	Guardian 2000
				
	Advanced Safety Features and Performance	All Around Versatility	Easy and Dependable	Entry-Level Simplicity
Functions	Heat and Stir	Heat, Stir, Heat and Stir	Heat and Stir	Heat, Stir, Heat and Stir
Plate Size	Round, 7" x 7", 10" x 10"	Round, 7" x 7", 10" x 10"	Round, 4" x 4", 7" x 7"	Round, 4" x 4", 7" x 7"
Display	5" LCD with Status Bar	5" LCD	3" LCD	Analog
Overheat Protection	●	●	●	●
SafetyHeat	●	●	●	
Temperature Probe	●	●	●	
Timer	●	●		
Data Transfer and Logging		●		
SmartRate	●	●		
SmartHeat	●	●		
SmartLink	●			
SmartPresence	●			
Programmability	●			
	Probes Suitability for Different Media			
	Stainless Steel Probes 30973654 (PT100, 25 cm) 30500590 (PT1000, 20 cm) 30500592 (PT1000, 25 cm)		Hastelloy Probe 30973655 (PT1000, 25 cm)	PTFE Probes 30500591 (PT1000, 20 cm) 30500593 (PT1000, 25 cm)
	Application Temperature Range			
	-40~400°C		-40~200°C	
	Aqueous Corrosion Resistance			
Organic Acids	A	A	A	
Alkaline & Salts	B	A	A	
Seawater	B	A	A	
Sulfuric Acid	B	A	A	
Hydrochloric Acid	D	A	A	
Hydrofluoric Acid	D	A	A	
Phosphoric Acid	C	A	A	
Nitric Acid	C	C	A	
Chlorine (wet)	D	A	A	
	A = Excellent B = Good C = Acceptable D = Not Recommended			

GUARDIAN SERIES

Accessories*




Vial and Test Tube Blocks



Base Plate and Handles



Temperature Probe

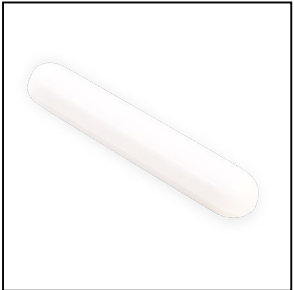

Spin Bars



Ultra Flex Support Kit



In Use Cover



Support Rod and Clamp Kit



Vessel Clamps

*Over 40 accessories available for Guardian Series



GUARDIAN 7000



GUARDIAN 5000



GUARDIAN 3000



GUARDIAN 2000

OHAUS Corporation

Headquartered in Parsippany, NJ, OHAUS Corporation manufactures an extensive line of balances and scales, lab equipment and lab instruments that meet the weighing, sample processing and measurement needs of multiple industries. We are a global leader in the laboratory, industrial and education markets, as well as a host of specialty markets, including the food preparation, pharmacy and jewelry industries. An ISO 9001:2015 manufacturer, OHAUS lab balances, industrial scales, lab equipment and lab instruments are precise, reliable and affordable, and backed by industry-leading customer support.

80776559_J 20240918 © Copyright OHAUS Corporation

OHAUS CORPORATION
7 Campus Drive
Suite 310
Parsippany, NJ 07054 USA

Tel: 800.672.7722
973.377.9000
Fax: 973.944.7177
www.ohaus.com

The management system governing the manufacture of this product is ISO 9001:2015 certified.



Ingeniously Practical