


Reveal® for Multi-Treenut

Rapid multiple allergen detection

Simple food allergen control

Reveal® for Multi-Treenut is a unique and innovative lateral flow test for rapid allergen detection of six tree nuts with one test. Reveal for Multi-Treenut is designed to screen environmental swabs and rinse waters for the presence of any one or a combination of almond, hazelnut, pecan, walnut, cashew and pistachio.

Our test detects allergen levels at low parts per million (ppm) and is perfect for on-site use with results in only 10 minutes.

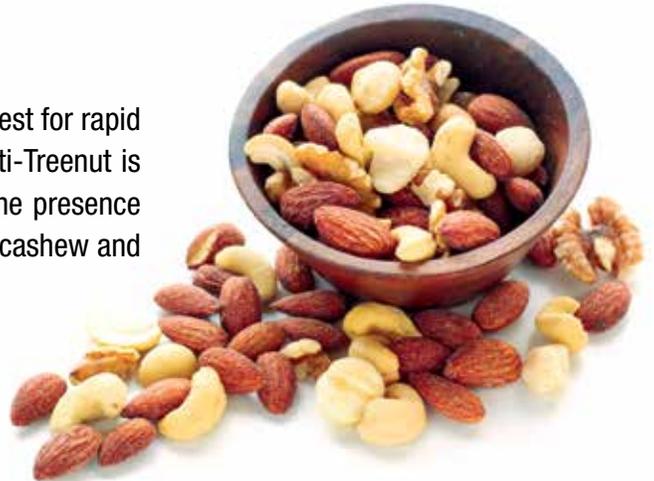


DETECTS

- ✓ Almond
- ✓ Walnut
- ✓ Pistachio
- ✓ Hazelnut
- ✓ Cashew
- ✓ Pecan



A new way to test

- Multiple allergens in one test → Cost and time effective
- Only test of its kind on the market → Unique
- Easy to use → No training required
- Rapid screening → Real-time decisions

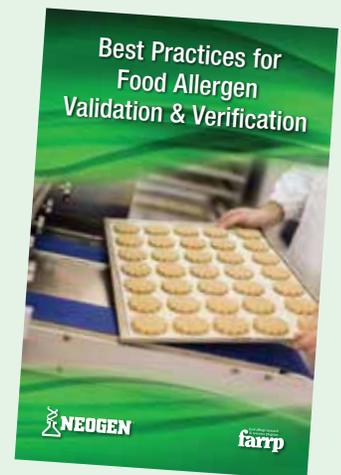

Contact us for a free Best Practices for Food Allergen Validation & Verification handbook



800-234-5333 (USA/Canada) or 517-372-9200
foodsafety@neogen.com • www.neogen.com



Reveal[®] for Multi-Treenut Procedure

Note: Please read kit instructions completely before performing test.

Rinse Sample Extraction



1. Open Type 11 extraction buffer sachet and add to a sample tube.



2. Add 1 mL of sample to the sample tube.




3. Secure cap and shake for 1 minute.

Swab Sample Extraction



1. Open Type 11 extraction buffer sachet and add to a sample tube.

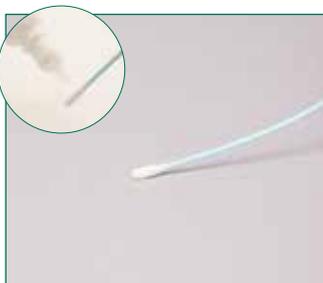

2. Gather the sample. *For dry surfaces moisten with swab wetting solution. Do not premoisten for wet surfaces.*



3. Return the swab to the extraction buffer and break off into the tube.



4. Secure cap and shake for 1 minute.

Test Procedure



1. Remove the lid and fill with liquid from the tube.



2. Dip the Reveal device into the liquid in the lid; ensure that the cavity is saturated with the liquid.



3. Leave the cavity saturated until the liquid is seen running into the test window.



4. Place the device on a flat surface and allow the test to develop for 10 minutes.



5. Interpret the results.

#8555

Reveal for Multi-Treenut

Rinse: 5-10 ppm

Swab: 5-20 µg/100 cm²
on Teflon

Testing time: 10 minutes

Materials Provided

10 lilac test devices

10 sample tubes and caps

10 sachets of Type 11 extraction buffer

10 sterile swabs

1 bottle swab wetting solution

Call **800/234-5333** to order or
visit us online at www.neogen.com



800-234-5333 (USA/Canada) or 517-372-9200
foodsafety@neogen.com | www.neogen.com

