Autoloop PRO

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Flame time
Precise settings of the required flame sterilization time

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Cooling period
The cooling time can be set for the inoculation loop material used - exact to the second

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Sterilization monitor
An intelligent sensor monitors the safe sterilization. Three setting options / off switch

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Easy to use

All parameters are easy to set; quick start button for immediate start

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Mount safely

Coordinated mounts and reliable optical detectors for inoculation loop holders

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Optimal sterilization

Sterilization monitor for safe sterilization of inoculation loops

Autoloop^{PRO}

The fully automatic carousel for flame sterilizing inoculation loops.

The inoculation loop carousel **Autoloop^{PRO}**, which is suitable for up to 4 inoculation loops, was developed to comply with current requirements in microbiological, cellbiological and biotechnological laboratories.

Modern design

The stable housing enables comfortable and easy access to inoculation loops. Removal positions on both sides make the carousel equally suitable for right and left handers.

More comfort

Keep all functions in view with the fully graphic display. The graphic menu control makes it easy to select all functions quickly. Flaming and cooling time can be adjusted to the second. During the flame sterilizing, the display shows the remaining flaming time. The **Autoloop^{PRO}** rotates and controls flaming automatically.

No unintentional use of hot inoculation loops: When flaming is completed, the display shows the remaining cooling time and the removal positions of cool inoculation loops. All the suspended inoculation loops can be flamed as needed in one full rotation of the carousel.

For additional flame sterilization (e.g. of bottle necks or test tubes) a foot pedal* can be activated during the automatic operation. For this purpose the **Autoloop^{PRO}** turns the inoculation loop out of the flame position and enables additional flame sterilization.

More safety

The nearly perpendicular position of the inoculation loop holder protects even the sleeve nut of the holder from possible contaminations. Additionally, an intelligent sensor of the **Autoloop^{PRO}** monitors safe sterilization.

Saves energy, time and costs

Precise flame sterilization time extends the life of the inoculation loops and reduces gas consumption. Continuous working during the flaming and the cooling phases of the inoculation loops saves a great deal of time and makes it possible to work efficiently.

Robust

Fabricated entirely of stainless steel, anodized aluminium and a display, protected by heat resistant glass, the **Autoloop^{PRO}** can withstand extreme laboratory conditions.

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Start options

Start immediately or sterilize all inoculation loops in a full rotation

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Everything visible at a glance

Removal positions of cool inoculation loops or the remaining cooling time, residual heat display and sterilization time

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Additional flame sterilization

of bottle necks or test tubes during the automatic operation by activating a foot pedal.

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The range:

Autoloop^{PRO}

includes 4 inoculation loop holders
4 inoculation loops
Pedestal
Foot pedal connector
Instruction manual
2-year warranty
Laboratory gas burner **not**
included

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Spray protector for Autoloop^{PRO}
for more safety when handling with
pathogenic material

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Spare inoculation loops
made of special stainless steel,
wire Ø 0.6 mm

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Technical data:

Technology:	Microprocessor
Display:	fully graphic dot-matrix
Rotational direction:	left / right, depends on removed holder
Inoculation loop holder:	1 - 4, with optical holder detection
Time of flame sterilizing:	1 - 15 sec
Time for cool-down period:	1 - 45 sec
Additional flame sterilization:	by foot pedal (optional)
Safety features:	temperature control flame activation warning indicator of remaining cooling time tilt sensor, residual heat display
Adjustable sterilization monitor:	1 - Off 2 - low intensity 3 - middle intensity 4 - high intensity
Stand by:	automatic unit switch off: 1 h (or with stand-by time of the burner)
Operating voltage:	via gas burner (foot pedal socket)
Casing:	stainless steel & aluminum, UV and solvent resistant
Measurements:	Pedestal (Art.-No. 8.000.400): 135 x 180 mm (w x d) Pedestal (Art.-No. 6.000.400): 135 x 135 mm (w x d) Height Autoloop ^{PRO} : 305 mm
Weight:	Art.-No. 8.000.400: 1200 g Art.-No. 6.000.400: 1100 g
CE:	EN 61326-1, EN 61010-1, EN 61010-2-010
EU guidelines:	2014/35/EU, 2014/30/EU, 2011/65/EU

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WLD-TEC GmbH